



# **PURELAB® Chorus 1 Complete\***





# **Purification Pack**

**ELC275** 

Lifetime expectancy\*\*: 6 months What is purified from the water?

Impact of not changing?

Weakly bound ions are released back into the water reducing resistivity



#### **ELC210** UV Lamp 185/254 nm

Lifetime expectancy\*\*: 12-18 months What is purified from the water?

Purifies TOC, inactivates DNA and RNA Impact of not changing?

Uncontrolled bacterial contamination and TOC levels are not reduced



#### Disinfection tablets

ECT1

Lifetime expectancy\*\*: 6 months What is purified from the water?

Sanitizes tubing or pipe work to remove bacteria, reduce bio film and protect the system against bacterial growth



### Pre-Treatment

**ELC241** 

Lifetime expectancy\*\*: 6 months What is purified from the water?

Chlorine and particulates

Impact of not changing? Chlorine, chloramine and organics remain increasing the burden on downstream consumables which will need more frequent and expensive changes



# Reverse Osmosis

**ELC240** 

Lifetime expectancy\*\*: 24 months What is purified from the water?

Majority of organics and inorganics Impact of not changing? High levels of organics

and inorganics remain in the water reducing the functional life of the DI Pack (Purification Pack) Note: Replacement need technician support



# Composite Vent Filter

**ELC216** 

Lifetime expectancy\*\*: 6 months What protection does it provide? Reduces CO<sub>2</sub> and airborne organics from entering the sealed system

Impact of not changing?

Increase in airborne contaminants and potential recontamination of purified water



# Point Of Use Filter\*\*\*

**ELC145** 

Lifetime expectancy\*\*: 3 months What is purified from the water? Particulates and bacteria Impact of not changing?

Potential recontamination of purified water and reduced flow rates

- \* A combination of treatments and technologies produce the final water specification
- \*\* Service life will depend on the application and feed water quality
- \*\*\* For DNA, RNA & endotoxin specification at point of use, please use a **ELC197 Biofilter**

